

AMBERLITE™ XAD™16HP N

Polymeric Adsorbent

Introduction AMBERLITE XAD16HP N is a polymeric adsorbent available in the form of white insoluble beads, supplied in the fully hydrated form. This resin has been prepared to meet high purity specifications. It is a non ionic, hydrophobic, cross-linked polystyrene copolymer, characterised by its unique macro reticular porosity, controlled pore size distribution and high surface area. AMBERLITE XAD16HP N has excellent physical resistance and thermal stability. In addition, it has a low swelling between solvent and aqueous media and can be used in column or batch operations.

Properties

Matrix	Macroporous cross-linked polymer
Physical form	White beads
Moisture holding capacity	62 to 70 %
Shipping weight	675 g/L
Specific gravity	1.015 to 1.025
Particle Size	
Harmonic mean size	0.600 - 0.750 mm
Uniformity coefficient	≤ 2.0
Fines content	< 0.300 mm : 3.0 % max
Coarse beads	> 1.18 mm : 5.0 % max.
Surface area	≥ 800 m ² /g
Porosity	≥ 0.6 ml/ml

Suggested Operating Conditions

Maximum operating temperature	150°C
Minimum bed depth	700 mm
Flow rate	5 to 40 BV*/h
Loading step usually	2 to 16 BV*/h
Washing step	1 to 2 BV/h
Regeneration step	1 to 2 BV/h
Rinse step	2 to 16 BV/h
Regenerants Methanol or other water miscible organic solvents (ethanol, acetone, Isopropanol, etc.)	
Dilute bases for weakly acidic compounds	
Dilute acids for weakly basic compounds	
Water where adsorption is from an ionic solution	
Hot water or steam for volatile materials	
Backwash Backwashing with water at 1 m/h and 15°C will produce a bed	
expansion of 70 %.	

* BV (Bed Volume) = 1 m^3 solution per m^3 resin

Food processing

AMBERLITE XAD16HP N meets the requirements of FDA 21 CFR 173.65 for removal of organic substances from aqueous foods except carbonated beverages. It can therefore be used for fruit juice decolourization and debittering and for the recovery of anthocyans, where the FDA regulation is applicable.



As governmental regulations concerning the use of resins for food processing applications may vary from country to country, it is recommended that potential users contact Rohm and Haas in order to assess the best choice of resin and optimum operating conditions.

In general, AMBERLITE XAD16HP N is ready to use after a simple regeneration followed by a rinse with 20 bed volumes of potable water.

Other applications In general AMBERLITE XAD16HP N is used for the adsorption of water soluble organic substances but it is not limited to this area.

AMBERLITE XAD16HP N is particularly useful in the separation and purification of antibiotics, vitamins, steroids, amino acids, enzymes, polypeptides, etc. in the pharmaceutical and organic chemical industries

For more information about DOW™ resins, call the Dow Water & Process Solutions business:

North America:	1-800-447-4369	
Latin America:	(+55) 11-5188-9222	
Europe:	+800-3-694-6367	
Italy:	+800-783-825	
South Africa:	+0800 99 5078	
Pacific:	+8007776 7776	
China:	+400 889-0789	
http://www.dowwaterandprocess.com		

Notice: No freedom from infringement of any patent owned by Dow or others is to be inferred. Because use conditions and applicable laws may differ from one location to another and may change with time, Customer is responsible for determining whether products and the information in this document are appropriate for Customer's use and for ensuring that Customer's workplace and disposal practices are in compliance with applicable laws and other governmental enactments. The product shown in this literature may not be available for sale and/or available in all geographies where Dow is represented. The claims made may not have been approved for use in all countries. Dow assumes no obligation or liability for the information in this document. References to "Dow" or the "Company" mean the Dow legal entity selling the products to Customer unless otherwise expressly noted. NO WARRANTIES ARE GIVEN; ALL IMPLIED WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE ARE EXPRESSLY EXCLUDED.

